

## Removal of Microspectrophotometer

1. Switch off PC and lamps
    - a. Note that lamp fan should be left on
    - b. Do not move DH-2000 until lamps are cool
  2. Disconnect fibre optics and pack PC
    - a. re-cap all ends
  3. Disconnect all cables and store
    - a. USB
    - b. RS232
    - c. Power cables
    - d. Ethernet
    - e. Security
  4. Pack HR2000 in flight case
  5. Remove cryo and goniohead
  6. Unmount microspec lenses
    - a. Remove springs and carefully extract lenses making sure not to drop bearings and pack
  7. Unmount stage
  8. Make sure packing foam is in position and lenses are protected
  9. Remount goniohead and reposition cryostream
  10. Store all bolts and tools
  11. Confirm that all parts are present as in Section 1
- Final check for cryostream position